

## Brief installation guide for bMCA software under Windows 8

BrightSpec, Niel, Belgium



03/2013

The installation of bMCA device (drivers) and software might fail under Windows 8 due to the new restricted policy for installing third party drivers (\*.INF) that does not contain digital signature.


To proceed with this driver installation you must disable the Windows 8 driver enforcement option.

The following steps show you how to do this:

Using the PC Settings, **press SHIFT key** while clicking at the Restart menu option from POWER

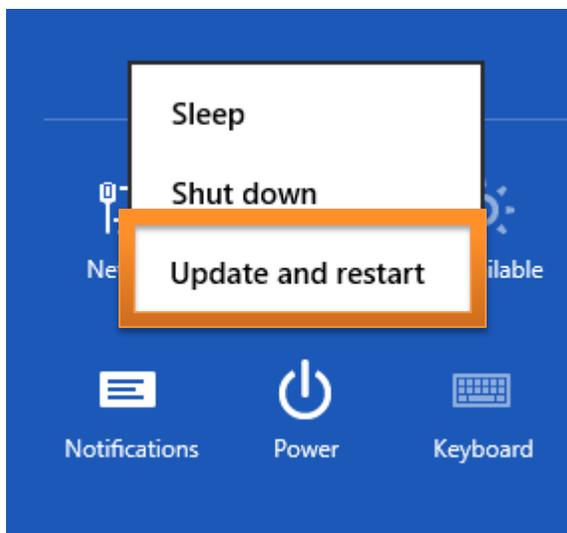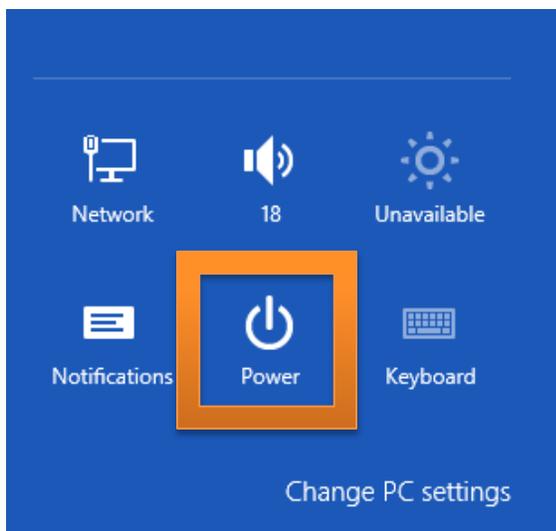

The following screen should appear

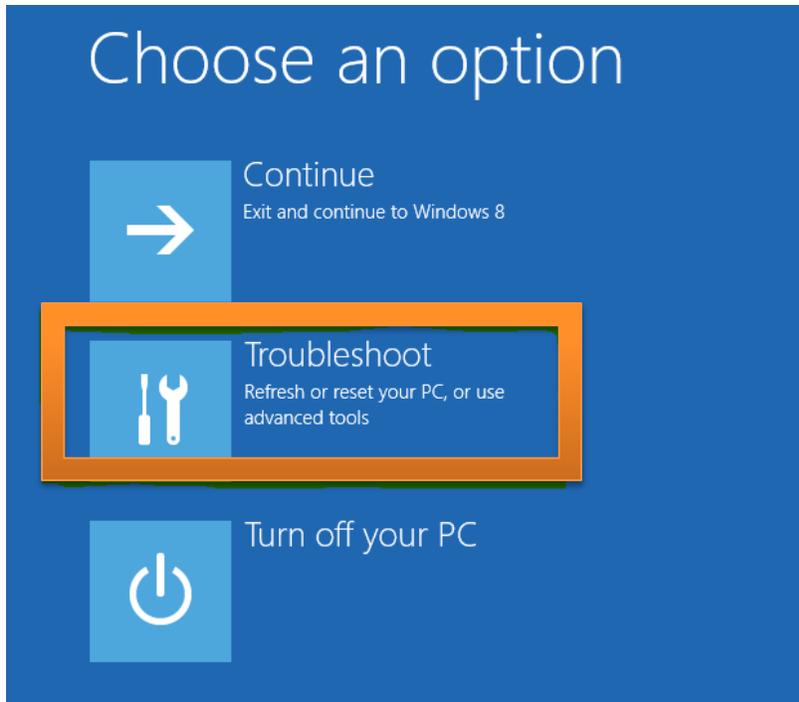


Chose the option “Troubleshoot”


From the Troubleshoot menu chose “Advanced options”, as shown below:



And from Advanced options select “Startup settings”

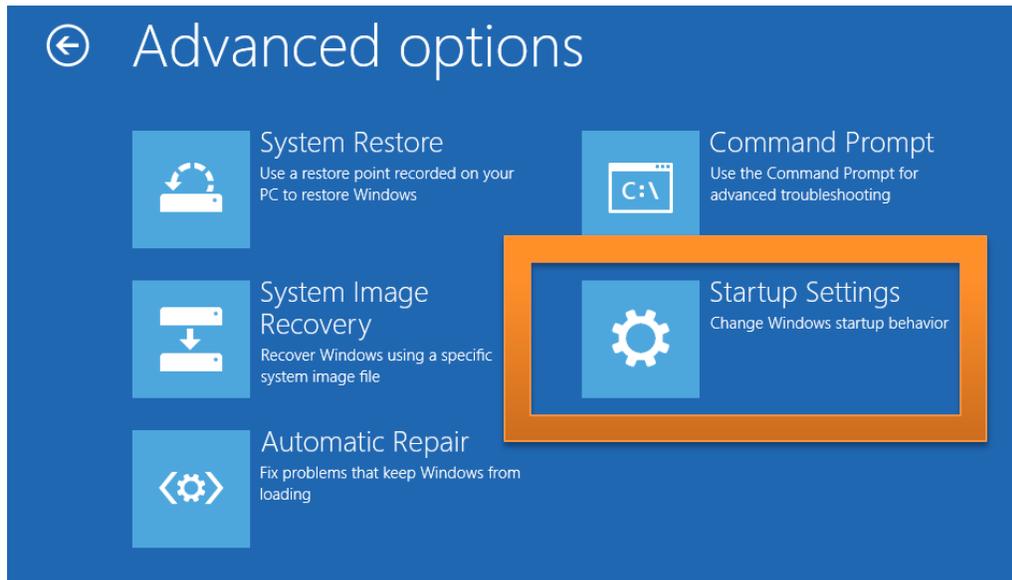

From “Startup Settings” screen chose Restart

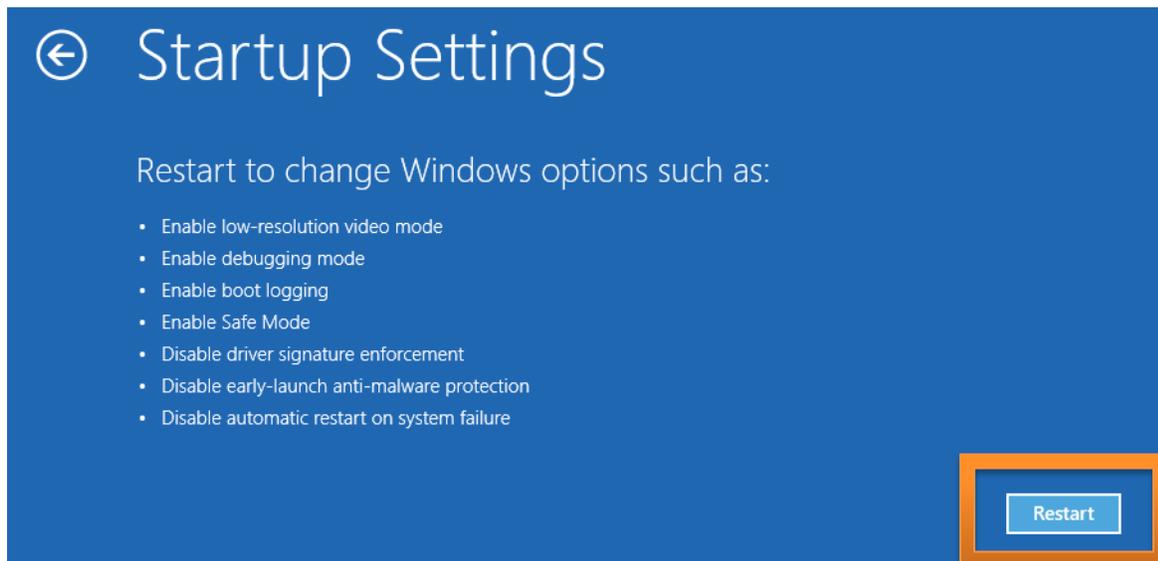

On next booting screen chose option 7 “Disable driver signature enforcement” by pressing “F7” key from the keyboard

# Startup Settings

Press a number to choose from the options below:

Use number keys or functions keys F1-F9.

- 1) Enable debugging
- 2) Enable boot logging
- 3) Enable low-resolution video
- 4) Enable Safe Mode
- 5) Enable Safe Mode with Networking
- 6) Enable Safe Mode with Command Prompt
- 7) Disable driver signature enforcement
- 8) Disable early launch anti-malware protection
- 9) Disable automatic restart after failure

Press F10 for more options

Press Enter to return to your operating system

After this the driver and software installation should progress as expected.